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Microscope Illuminator Hal
with Reflector

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Microscope Illuminator Hal with Reflector

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
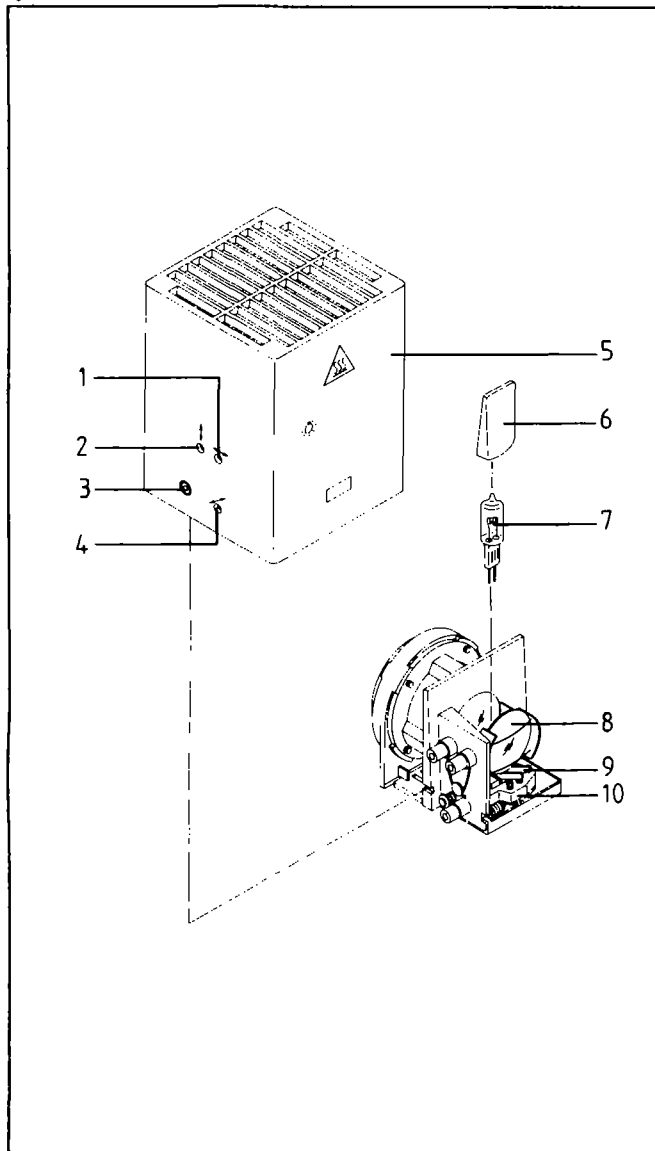
The microscope illuminator Hal (44 72 17) consists of the lamp housing Hal with collector, reflector, adjustable lamp socket, heat protection filter and the halogen lamp.

Replacement and adjustment of the halogen lamp are described in this instruction manual. Attachment and connection to the microscopes Axioplan, Axiophot, Axiotron, Axioskop and Axiovert are included in the microscope instructions.

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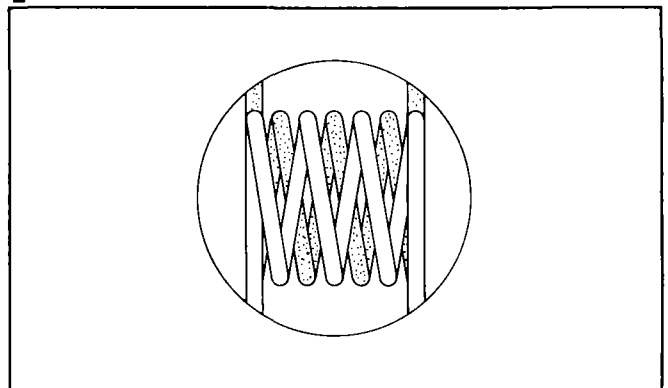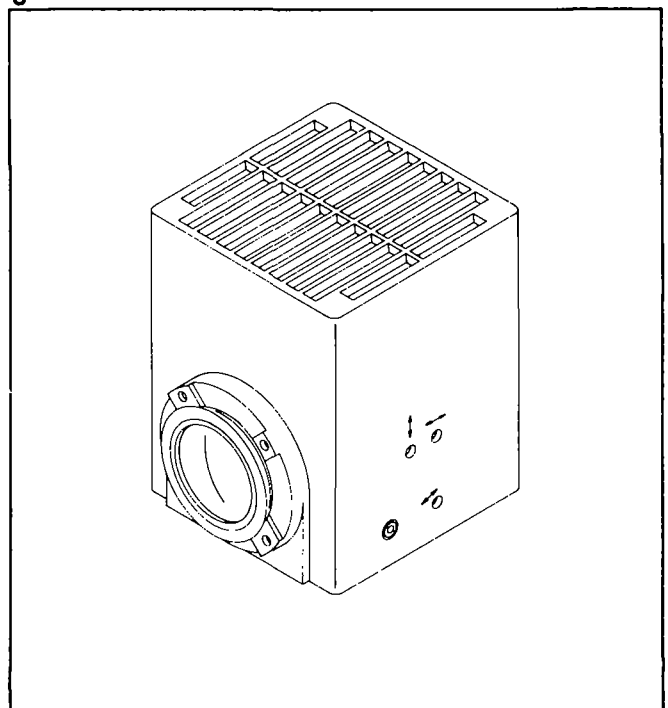
Replacement of the halogen lamps 12V 100W (380079-9540) and 12V 50W (380079-9550)

This warning sign at the back of the illuminator means: **Caution: hot surface!** Do not touch until cool.

- Switch off microscope before lamp replacement and pull out plug from lamp supply unit. Wait until lamp is cool. The illuminator can remain mounted on the microscope for the lamp exchange.
 - Loosen screw (3) and lift off lamp housing (5).
 - Pull defective lamp bulb out of lamp socket (10) by simultaneously pressing down both springs (9).
 - Insert new lamp (7) using protective sleeve (6) as follows:
 - Pull lamp out of protective sleeve until pins are fully visible (avoid fingerprints).
 - Hold top of lamp using protective sleeve and insert lamp into the socket while the springs (9) are pressed down. Press down springs again to enable automatic positioning of the halogen lamp. Attach lamp housing again and fix it using screw (3). Connect illuminator cable to power supply socket again and switch on the lamp (see operating instructions of the microscope).
- Fix adhesive label supplied with the halogen lamp to the frame at the back of the lamp housing.

Adjustment of the halogen lamp with reflector

- The Hal microscope illuminator is attached at the microscope.
- Remove diffusion disk and filter, if used, from light path (see microscope instructions).
- Focus on specimen using an objective with magnification 40x or higher. Find a vacant specimen position. Fully open aperture diaphragm in condenser. When the eyepiece is removed and the centering telescope inserted into the tube, view coil image and reflector image and check whether the gaps in the reflector image are covered (see Fig. 2). If not, make correction as follows:
 - Image the lamp coil sharply using screw (1), use screw (2) for vertical adjustment and screw (4) for horizontal adjustment. Adjustment is correct when the gaps in the reflector image are fully covered. The fixed reflector (8) remains permanently adjusted.
 - Insert eyepieces into the tube and replace diffusion disk and filter, if used, into the light path again.

Check homogeneous illumination of pupil image and correct it, if required, using screw (1).

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Notes on safety and maintenance

Always switch off the microscope and pull out the lamp plug before you open the illuminator and wait until lamp is cool.

The halogen lamp must always be operated in the closed lamp housing to avoid burning the hands and dazzling the eyes.

Changes and/or repairs of the instrument should be carried out only by the manufacturer or persons authorized by him to do so.

Use only the halogen lamps mentioned above. Other commercial lamps will cause optical and perhaps also mechanical difficulties.

Manufacture, inspection, assembly, maintenance and repair are made in compliance with German and international instructions (Good Manufacturing Practice).

Clean painted surfaces using a cloth soaked in a mixture of water and detergent and dry with a lint-free cloth. Do not use solvents.

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